

Installation Guide

Visagraph® III Standalone

Updated April 18, 2008



REQUIREMENTS


Please check the following table to ensure that you meet the minimum system requirements before continuing on to installation.

	Visagraph® Standalone Workstation Requirements
OS	Win 2000 Pro, XP
CPU	Pentium 1.0+ Ghz
RAM:	512 MB (750 MB recommended)
Hard Disk Space	25 MB free + 50Kb per recording
Interface	USB Port
Monitor Resolution	1024 X 768
Browser	IE 6.0+ w/ Java Sun Plug-in 1.5.0_5+

INSTALLATION INSTRUCTIONS


Windows 2000/XP/Vista

1. Place the CD Into the CD-ROM Drive and wait for the CD to start. If the CD doesn't start, open the CD Drive and double click the Visagraph.exe file.
2. At the *Welcome to the Visagraph Setup Wizard* screen, click the *Next* button.

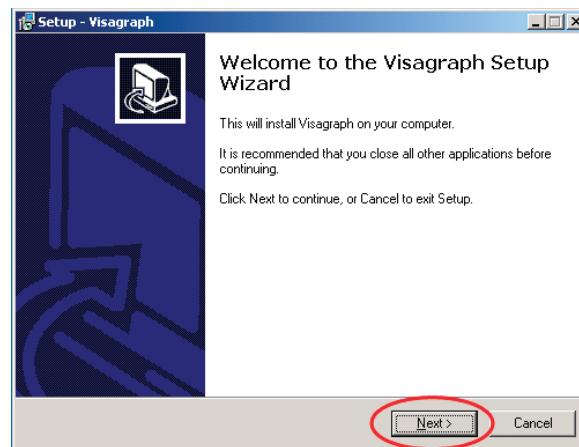

3. Select the port you would like the Visagraph® III to use. (The default port is 80. If you would like to use another port for your Visagraph® III, enter it into the field and click *Check Availability*.) If the port is not in use, click *OK*.

5. Choose your destination path and click *Next*.

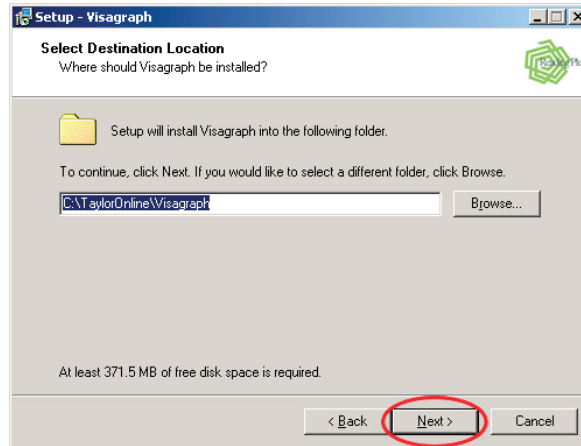

6. If you are running Windows XP sp2, you might be presented with a screen asking to accept Apache. Click *Unblock*.



Once the installer is complete, click the *Finish* button. You may now begin using the Visagraph® III.